

Food and Environmental Proficiency Testing Unit

The final date for return of results is indicated on the accompanying request/report form

Contact details:

The Organisers - FEPTU
UK Health Security Agency
61 Colindale Avenue,
London, NW9 5EQ, UK.
Tel: +44 (0) 20 8327 7119
e-mail: foodega@ukhsa.gov.uk



Results to be returned via an online survey as per email sent to your laboratory

Mycobacterium spp. - Food and Environmental Proficiency Testing Unit

Mycobacterium spp. in Water Scheme – Instruction Sheet

The LENTICULE® discs are provided in screw-cap plastic vials (with a desiccant). Each LENTICULE disc rests on the filter support containing the desiccant. The **desiccant** should be orange in colour. Please contact us if this is not the case. Each LENTICULE disc requires reconstitution by a process of rehydration and dispersion prior to examination, as described below.

Refer also to the Safety Data Sheet: Safety Data Sheet - LENTICULE

Storage:

- Store the samples at -20 ± 5°C on receipt.
- Allow the LENTICULE discs to reach ambient temperature (5 10 minutes) before reconstituting in diluent.

Reconstitution:

- Open the sample container and transfer the LENTICULE disc into approximately **9mL** 0.1% peptone saline (maximum recovery diluent (MRD)) that has been allowed to reach ambient temperature.
- Leave to stand at ambient temperature for 10 -12 minutes to rehydrate.
- Tighten the cap of the bottle and shake to disperse the micro-organisms.
- Transfer all the inoculated MRD to a sample bottle containing 1L sterile deionised or distilled water at ambient temperature.
- Rinse with approximately 2mL from the sample bottle, ensuring all liquid is transferred back to the 1L sample.
- Disperse the inoculum by inverting approximately 30 times.

Examination:

- Each reconstituted sample is equivalent to 1L water.
- Undertake the sample examinations between 30 45 minutes of reconstitution.
- Examine in accordance with your routine procedures.

Return of Results:

- Return your results on-line using the web-based reporting system.
- The distribution will be closed on the deadline date for return of results.
- Results cannot be accepted after the distribution has closed.

FEPTU840.04 Page 1 of 1